**CENTRÍFUGAS  
CENTRIFUGES  
CENTRIFUGEURS**

**Analógicas / Analogical / Analogiques**

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Thank you for choosing this equipment. We sincerely wish that you enjoy your analogical Nahita centrifuges. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:

- ◆ This manual is inseparable from the analogical centrifuges Nahita, so it should be available for all the users of this equipment.
- ◆ Follow all required applicable laws about safety and health (electrical security devices, inflammable, toxic and/or pathological products, etc.).
- ◆ You should carefully handle the analogical centrifuges avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the analogical centrifuges to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the analogical centrifuges.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to (asistencia@auxilab.es).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the analogical centrifuges warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the analogical centrifuges warranty.
- ◆ Neither fuses (1A) nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

**INDEX OF LANGUAGES**

Spanish .....	2-8
English .....	9-15
French.....	16-22

**INDEX OF CONTENTS**

1. USES OF THE INSTRUMENT .....	10
2. BASIC THEORY.....	10
3. DESCRIPTION .....	12
4. TECHNICAL SPECIFICATIONS .....	12
5. INSTALLATION / SETTING UP.....	11
6. MAINTENANCE AND CLEANING.....	14
ANNEX I: CE CERTIFICATE .....	23

**1. USES OF THE INSTRUMENT**

Compact centrifuges, they sturdy in design and can be used for a variety of applications.

Models with mechanical regulation have head for tubes of different volumes and with system of elastic suspension.

There are some examples of different activities in which this centrifuge is used as a technique of separation:

- a) Separation of serum or plasma.
- b) Cell concentration.
- c) Separation of substances.
- d) Biochemical laboratories.

**2. BASIC THEORYENGLISH**

The centrifuge is an instrument that provides a technique of separation that is based on the movement of the particles. In this way the particles are moved towards the end of the rotating axis depending on their different masses and shapes.

The Relative Centrifuge Force (R.C.F.) is the required force so that the separation can be carried out. The units of this force are expressed in number of times higher than the gravity (g).

- have a balance to adjust the weights of the different tubes.
- ◆ This is absolutely necessary since if the rotor is not well balanced it can produce shakings, if these shakings are small it is difficult to notice them, then little by little the centrifuge wears away and then the breaking of the tube holders increases and it produces a worse sedimentation. It is advisable to change all the tube holders at the same time.
  - ◆ Do not stop the centrifuge by the application of any external force, in this way you will avoid the possible mixture of the separate parts.
  - ◆ In case of damage or the change of the brushes, you have to get in touch with your supplier for the repair by the Nahita Technical Service.

***Cleaning***

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the analogic centrifuges and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.
- ◆ With high g-values, there is a possibility that tube breakage will occur. Should this happen, the centrifuge rotor, adapters and the rotor chamber must be thoroughly cleaned and all broken particles removed immediately.

**INSTRUCTIONS ON ENVIRONMENT PROTECTION**

At the end of its life cycle, please do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

correct use and intended purpose of the machine.

- ◆ You should put the Schuko standard wire in a horizontal plane stable table, having a safety area of at least 30 cm per side.
- ◆ Follow all required applicable laws about safety and health (electrical security devices, inflammable, toxic and/or pathological products, etc.).
- ◆ Do not place the analogic centrifuges near any warm supply (burners, blowlamps, etc), nor expose it directly to the sun. Avoid vibrations, dust and dry environments.
- ◆ When you are not using the analogic centrifuges for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the analogic centrifuges by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

## 6. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

*Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.*

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The analogic centrifuge is supplied with a Schuko standard wire and it should be connected to a current wire provided with an earth wire, it should be handy to be disconnected in case of emergency.
- ◆ The analogic centrifuge is supplied with a Schuko standard wire and it should be connected to a current wire provided with an earth wire, it should be handy to be disconnected in case of emergency.
- ◆ Never try to repair the analogic centrifuges by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the analogic centrifuges (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.
- ◆ Use tubes that are resistant to the R.C.F.
- ◆ Balance the centrifuge placing tubes that are equal in weight, shape and size in symmetrical positions. If it is necessary, we will use tubes filled with water when one of the holes is empty. In a centrifuge it is very important the concept of weight, then it is advisable to


The R.C.F. is calculated by means of the following formula:

$$F.C.R. = 1,118 * 10^{-5} * r * n^2$$

Being  $1,118 * 10^{-5}$  a constant.

r = horizontal distance of the radius (in centimetres) from the centre of rotation up to the bottom of the tube.


n = speed of rotation expressed in revolutions per minutes (r.p.m.).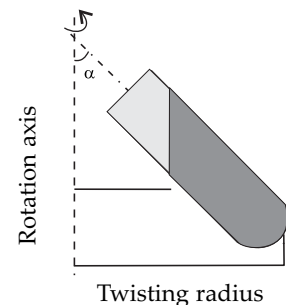


## 3. DESCRIPTION

- |                          |                           |
|--------------------------|---------------------------|
| 1. On/off switch         | 4. On/off indicator light |
| 2. Speed adjustment knob | 5. Lid                    |
| 3. Time adjustment knob  | 6. Rotor                  |


[Model 2615]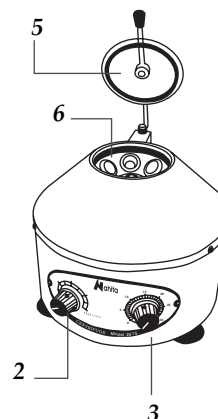

[Model 2690 and 2690/5]

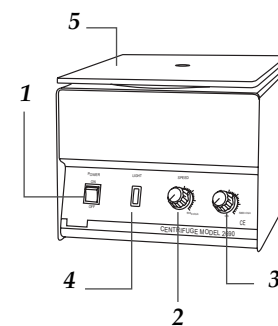[Model 2650]
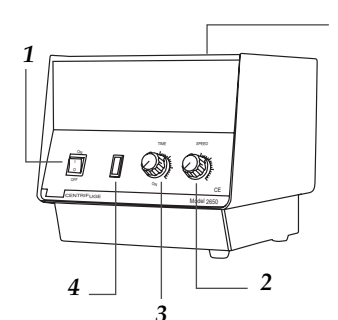
#### 4. TECHNICAL SPECIFICATIONS

Model	2615	2650	2690	2690/5
Control	Analogical			
Head	Inclined 45°			
Max. speed	4000 r.p.m	4000 r.p.m	5000 r.p.m	5000 rpm
Max. R.C.F	1790xg	2147xg	3130xg	3130xg
Capacity	6x15mL	12x15mL	8x15mL	16x5mL
Timer	0-30min			
Motor	Elastic suspension			
Power supply	220V-230V 50Hz			

#### 5. INSTALLATION / SETTING UP

##### Preliminary inspection

Unwrap the balance, take off the involving plastic and take off the poliespan protection in which it comes fitted.

Please make sure that it does not present any damage because of the shipment. If it does please communicate to your transport agent or suministrator immediately, so that they can make the claims in the correct time limit.

*We will only accept any equipment return within 15 days after delivery and provided it comes in its original package and with all accessories and documents supplied*

Please check that all the accessories are enclosed with the equipment:

- Tube holders
- Schuko standard wire
- Instruction manual
- Warranty

##### Installation

◆ Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

**PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT**

◆ Please put the analogical centrifuges on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. Do not put the analogical centrifuges near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.

- ◆ Avoid inflammable or toxic substances in the working area.
- ◆ The analogical centrifuges is supplied with a Schuko standard wire
- ◆ Please insert the wire that feeds the AC electric current in the base of current 220V 50Hz  $\pm 10\%$  provided with earth wire and to the other end to the analogical centrifuges connector.

*Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220V 50Hz  $\pm 10\%$ .*


If you are not using the analogical centrifuge for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

##### Starting the centrifuge

Open the cover:

- ◆ (According model) press the button which is in Right-lateral and open the cover Put the test sample into centrifuge tubes carefully.

**IMPORTANT: Make sure that the samples placed in the symmetrical tubes are of equal weight to avoid serious shaking due to big difference in weight, close the cover.**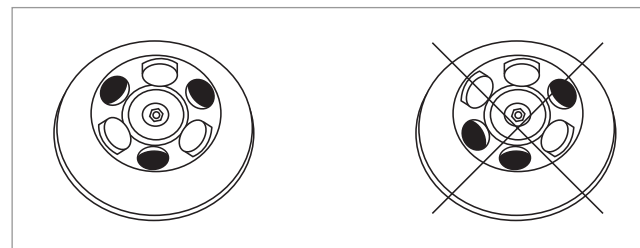

Power:

- ◆ Turn on the switch (1) and the light is lit to indicate that the power is connected. (Only on Model 2650, 2690, 2690/5)
- ◆ Adjust the speed and time with knob respective: speed knob (2) and Time knob (3).

Starting the centrifuge:

- ◆ When adjust the time to start the centrifuge. If you want to continuous run turn the knob in opposite clockwise to the stop limit (according model)

Stop

- ◆ Once it past times the rotor stop. If you want to interrupt a centrifuge run, turn the knob timer (3) in klokwise.

##### Security

- ◆ The centrifuge must not be operated by unqualified persons not familiar with the